

SensoScientific MAC ID & Probe SRS Numbers v.6/2026

How to locate the MAC ID on your OTA node and the SRS Number on your probe.

The OTA node is the model of SensoScientific data logger that displays the temperature and the Probe is what is in the unit monitoring temperatures and replaced annually. Probes are required to be certified and calibrated, certificates of calibration must be available if requested.

OTA Node

The MAC ID is located on the Node Display. The MAC ID consists of 12 numbers and letters. Each MAC ID is unique, each device has its own unique ID.

(For example: 509CB1779C30)

You may be asked for the MAC ID during troubleshooting and will need to provide it during annual re-enrollment.




SRS Probe:

The SRS number is located on two places on the cord of the probe.

The white tag will reflect the SRS Number only. The blue tag lists the SRS Number, Calibration Date and Calibration Due date.

Please note: the Calibration Due Date is one year from the calibration date. Certificates of Calibration can be located and printed/PDF saved on the [SensoScientific Cloud portal](#).

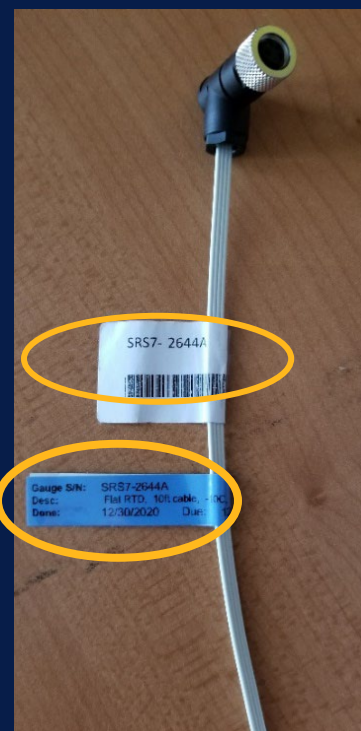

IMPORTANT: Do NOT contact SensoScientific directly. Facilities must direct all SensoScientific related questions to the Utah Department Health and Human Services, Immunization Program at vacteam@utah.gov.

This includes, but not limited to, requesting probes, adding and removing users and troubleshooting connection issues.